



# Fluorescence Filter Cubes

Images courtesy of : (1) Michael W. Davidson, National High Magnetic Field Laboratory, USA (2) Dylan Burnette, Paul Forscher Laboratory, Yale University, USA (3) Takeharu Nagai, Ph.D., Professor, Laboratory for Nanosystems Physiology, Research Institute for Electronic Science, Hokkaido University

Qdot, and the Q logo are trademarks of Quantum Dot Corporation.  
Texas Red® is a trademark of Molecular Probes, Inc. Cy® is a trademark of Amersham Biosciences, Ltd.  
Company names and product names appearing in this brochure are their registered trademarks and trademarks.

Monitor images are simulated.  
Specifications and equipment are subject to change without any notice or obligation on the part of the manufacturer. March 2010 ©2005/2007/2010 NIKON CORPORATION



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
The products detailed in the brochure are controlled by the Japanese Foreign Exchange and Foreign Trade Law and the International Export Control Regime. If there is a possibility that they may be utilized for the development of weapons of mass destruction, etc., they shall not be exported without authorization from the government.

**WARNING** TO ENSURE CORRECT USAGE, READ THE CORRESPONDING MANUALS CAREFULLY BEFORE USING THE EQUIPMENT.

**NIKON CORPORATION**  
Shin-Yurakucho Bldg., 12-1, Yurakucho 1-chome  
Chiyoda-ku, Tokyo 100-8331 Japan\*  
phone:+81-3-3216-2375 fax:+81-3-3216-2385  
<http://www.nikon.com/instruments/>  
\*from May 6, 2010

**NIKON INSTRUMENTS INC.**  
1300 Walt Whitman Road, Melville, N.Y. 11747-3064, U.S.A.  
phone: +1-631-547-8500; +1-800-52-NIKON (within the U.S.A. only)  
fax: +1-631-547-0306  
<http://www.nikoninstruments.com/>  
**NIKON INSTRUMENTS EUROPE B.V.**  
Laan van Kronenburg 2, 1183 AS, Amstelveen, The Netherlands  
phone: +31-20-44-96-300 fax: +31-20-44-96-298  
<http://www.nikoninstruments.eu/>  
**NIKON INSTRUMENTS (SHANGHAI) CO., LTD.**  
CHINA phone: +86-21-5836-0050 fax: +86-21-5836-0030  
(Beijing branch) phone: +86-10-5869-2255 fax: +86-10-5869-2277  
(Guangzhou branch) phone: +86-20-3882-0552 fax: +86-20-3882-0580

**NIKON SINGAPORE PTE LTD.**  
SINGAPORE phone: +65-6559-3618 fax: +65-6559-3668  
**NIKON MALAYSIA SDN. BHD.**  
MALAYSIA phone: +60-3-7809-3688 fax: +60-3-7809-3633  
**NIKON INSTRUMENTS KOREA CO., LTD.**  
KOREA phone: +82-2-2186-8410 fax: +82-2-555-4415  
**NIKON CANADA INC.**  
CANADA phone: +1-905-602-9676 fax: +1-905-602-9953  
**NIKON FRANCE S.A.S.**  
FRANCE phone: +33-1-4516-45-16 fax: +33-1-4516-45-55  
**NIKON GMBH**  
GERMANY phone: +49-211-941-42-20 fax: +49-211-941-43-22  
**NIKON INSTRUMENTS S.p.A.**  
ITALY phone: +39-055-300-96-01 fax: +39-055-30-09-93  
**NIKON AG**  
SWITZERLAND phone: +41-43-277-28-67 fax: +41-43-277-28-61

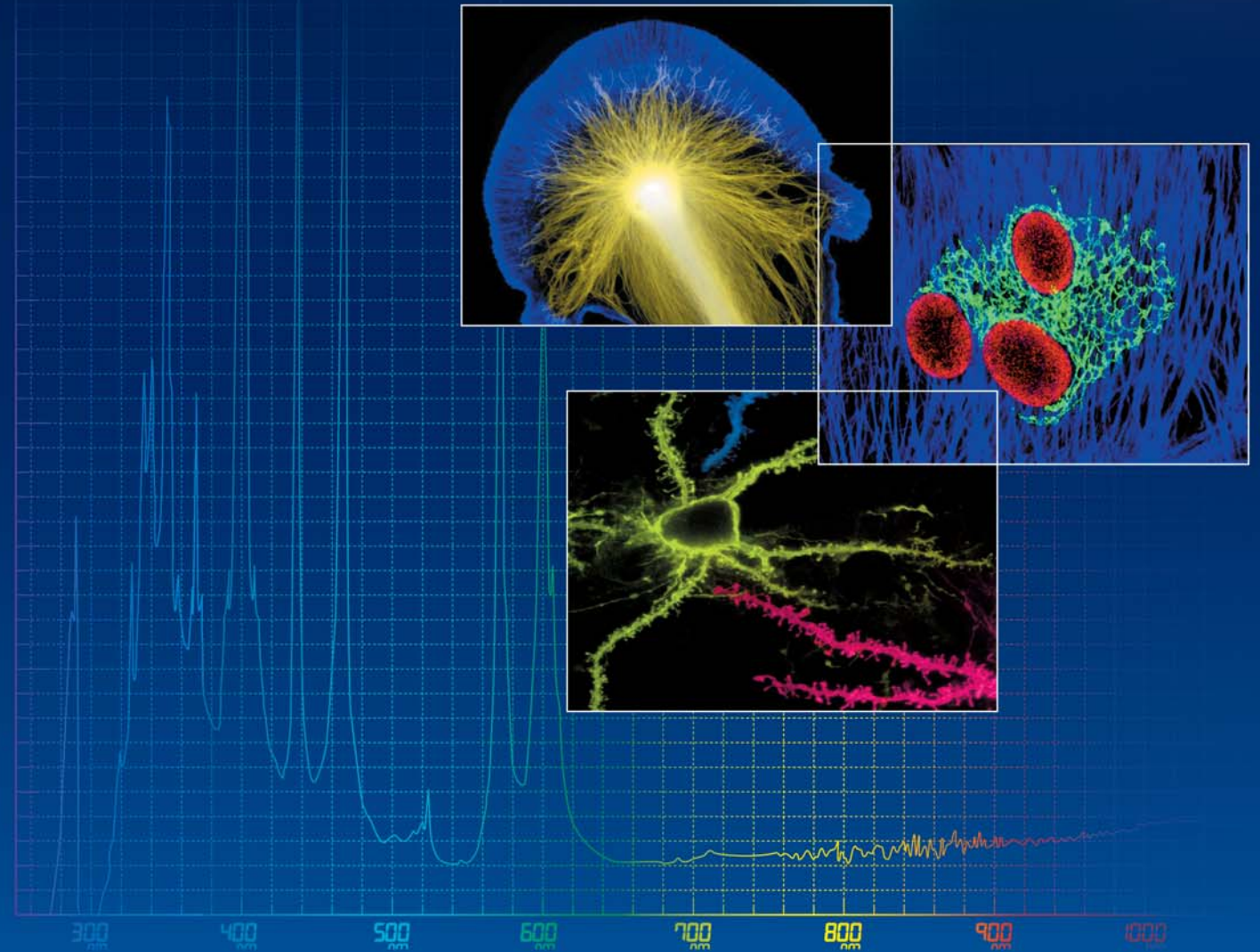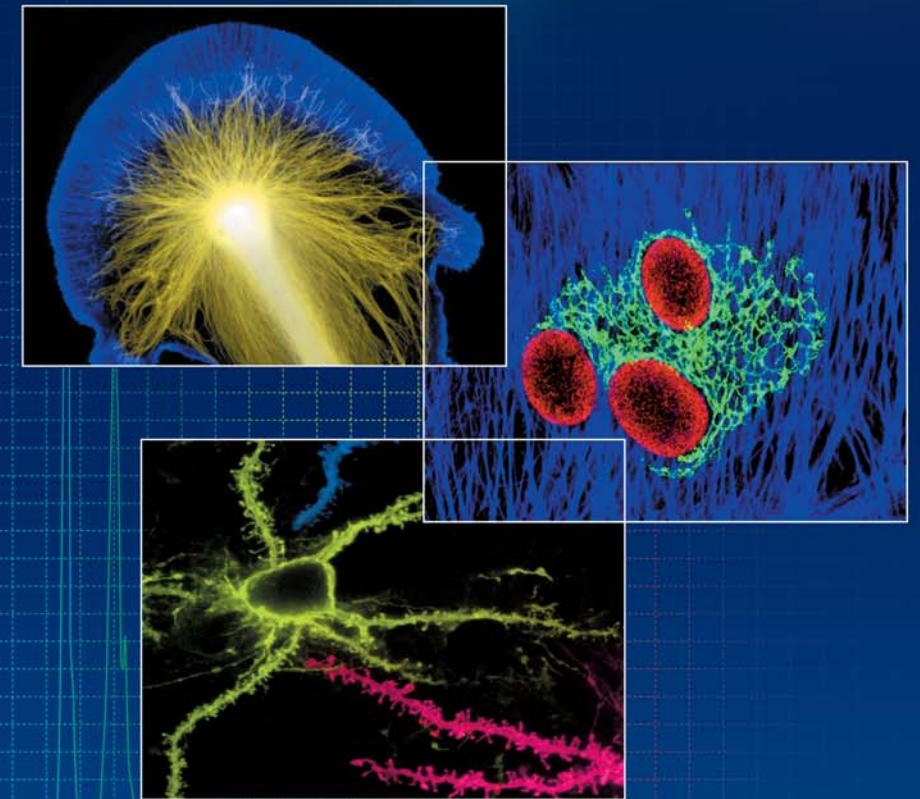



# Fluorescence Filter Cubes from Nikon




Nikon offers a wide range of filter cubes, from general epi-fluorescence use to those dedicated for a specific fluorescent reagent, to supporting today's variety of fluorochromes. With a standard filter diameter of 25 mm, commercially available fluorescence filters can be exchanged by users according to the desired use.

## Product Lineup

### Fluorescence filter cubes

General fluorescence filters corresponding to various excitation colors such as B-2A and G-2A. Since many of the barrier filters are the long-pass type, numerous reagents can be supported by a single filter.

#### Understanding filter names

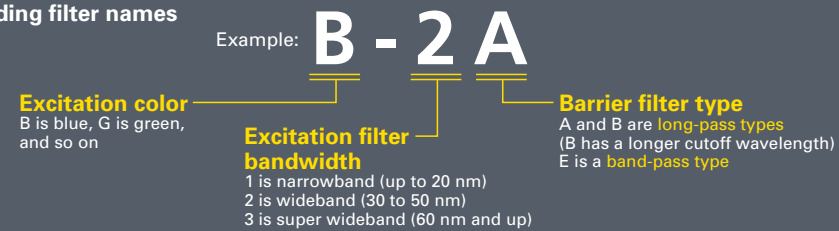

### Filter cubes for fluorescent reagents and fluorescent proteins


Filters corresponding to specific fluorescent reagents such as DAPI and FITC. Since the band-pass type is common on the barrier filter side, autofluorescence of plants, for example, is suppressed, enabling clear images with low background noise.

### High-quality filter cubes for fluorescent protein

The Wavelength cut-on and cut-off rise to peak is very steep, much more than for ordinary filter cubes for fluorescent proteins, thereby enhancing transmittance. Employing this filter makes it possible to obtain extremely clear, bright and non-overlapping fluorescent images.

### Multiband filter cubes


Filters that enable the simultaneous observation of double or triple staining techniques such as DAPI-FITC-Texas Red. Since there is no need to switch filter cubes in multi-stained fluorescence observation, there will be no position deviation of fluorescence filters, and no need to use the merge function of capture software when shooting with a color camera.

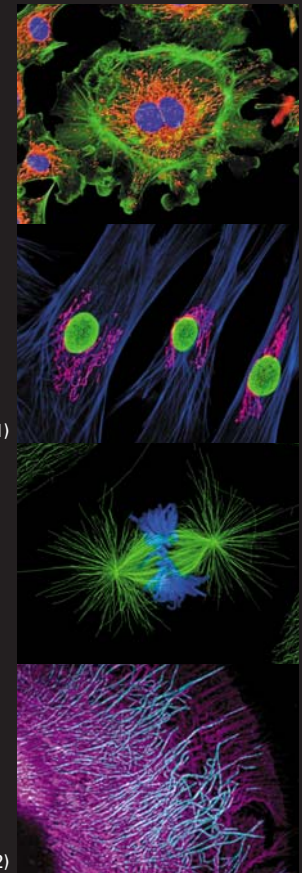

## Nikon Technology Maximizes the Potential of Fluorescence Images

### Noise Terminator

Nikon employs the proprietary Noise Terminator in its fluorescence systems to obtain even brighter fluorescence images. The Noise Terminator prevents excitation light (stray light) that barely passes through without being totally reflected by the dichroic mirror from scattering within the filter cube and having an effect on the image being observed. This makes it possible to obtain fluorescence images with extremely low background noise and a high S/N ratio.

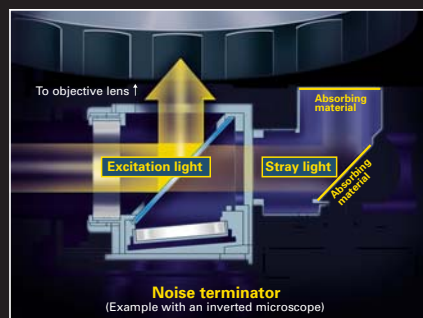

### High-performance filter cassette holder

The epifluorescence filter cassette holder for the Ti-E/-U/-S minimizes the deviation of the focal point of the light source on the objective pupil plane due to filter cassette switching. It is the optimum choice for TIRF illumination in all types of excitation methods.



# High Quality Filter Cubes GFP-HQ/CFP-HQ/YFP-HQ NEW




Enjoy bright, accurate detection, even for the samples with weak fluorescence. Nikon offers super high-performance/high-quality fluorescent filters for advanced research that demands precise signal detection.

### Main features of the HQ series of high-precision fluorescence filters


- Delivers wavelength characteristics with an extremely sharp rising edge.
- Delivers 93% to 97% transmittance (both excitation and barrier filters).
- Cuts excitation time nearly in half by increasing brightness by 50% or more.




## CFP-HQ *Caf1-SECFP-KDEL expressed in an endoplasmic reticulum*

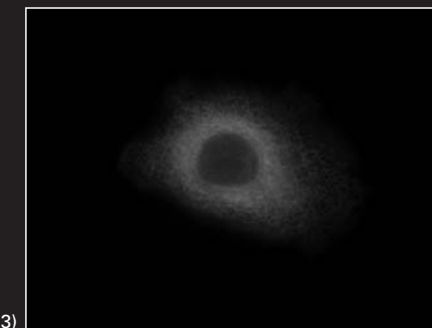

Using a normal filter

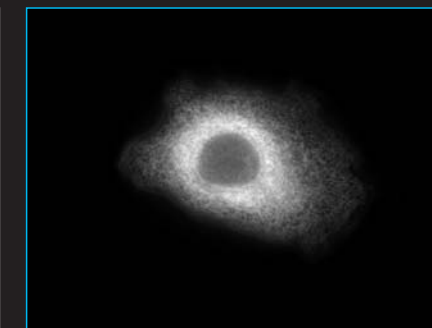


Using an HQ filter

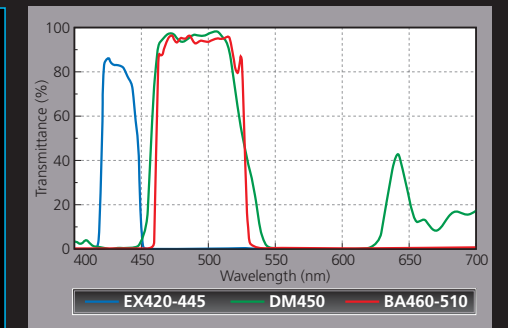


## GFP-HQ *SEGFP expressed in cytoplasm and nucleus*




Using a normal filter

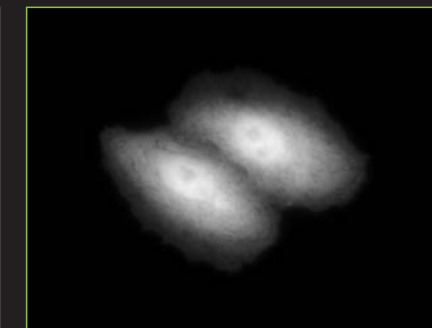


Using an HQ filter

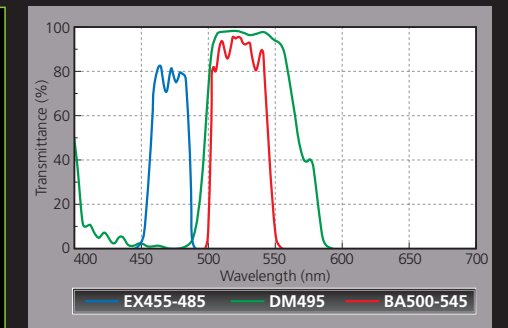


## YFP-HQ *CoxIV-Venus expressed in mitochondria*

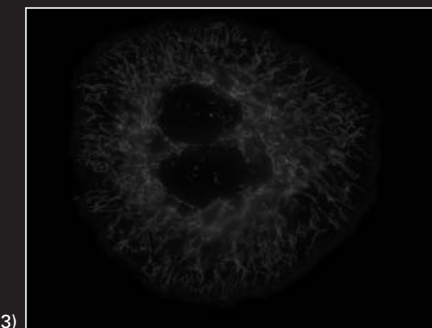

Using a normal filter

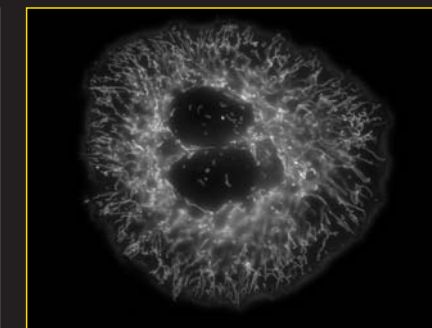

Using an HQ filter
















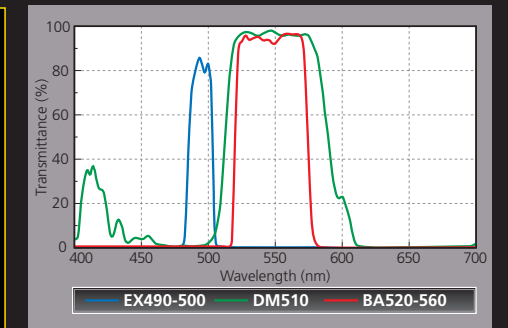


EX: Excitation filter DM: Dichroic mirror BA: Barrier filter

# Fluorescence Filter Cubes from Nikon



## Spectral characteristics table for filters

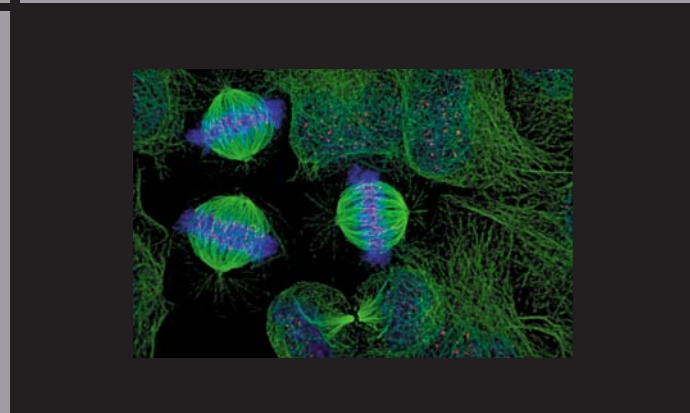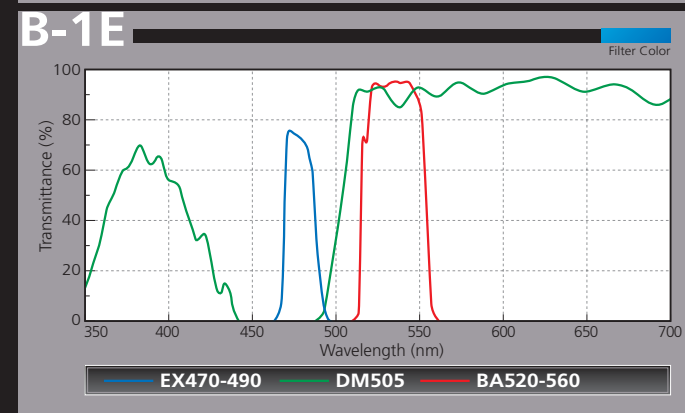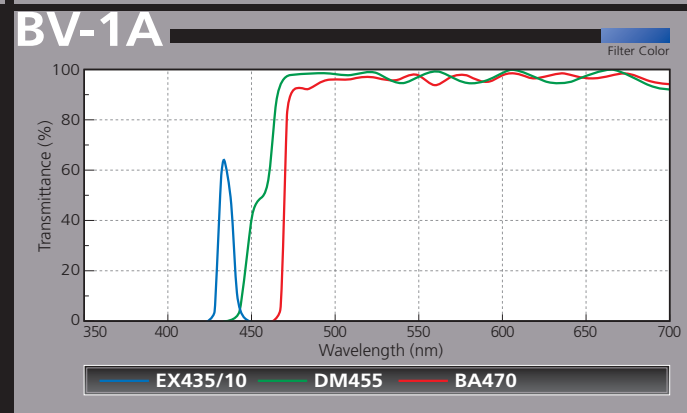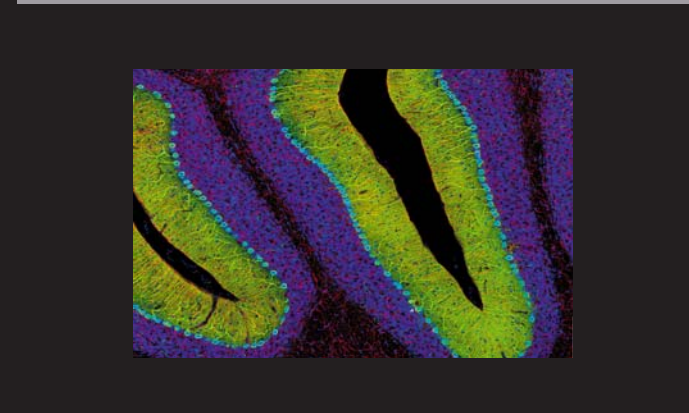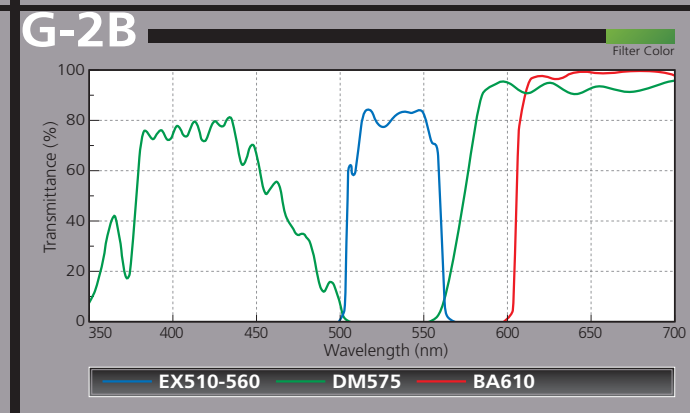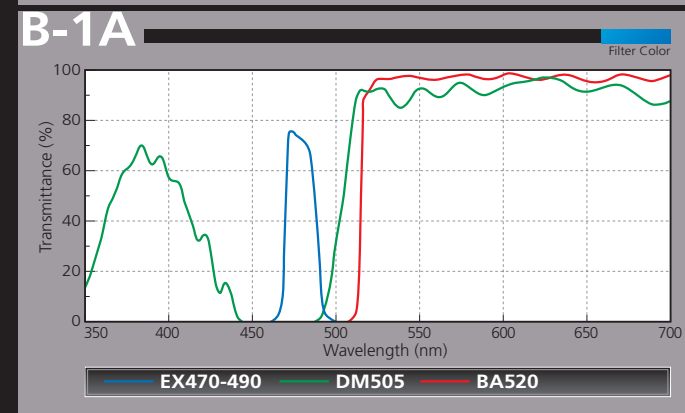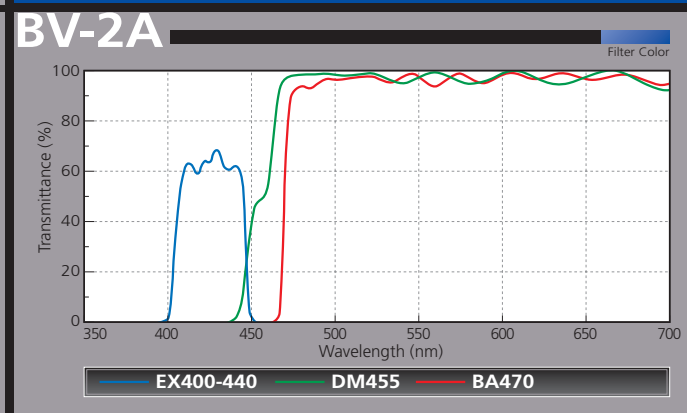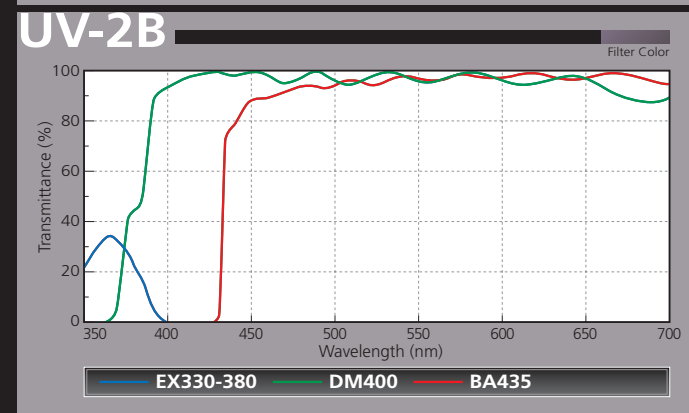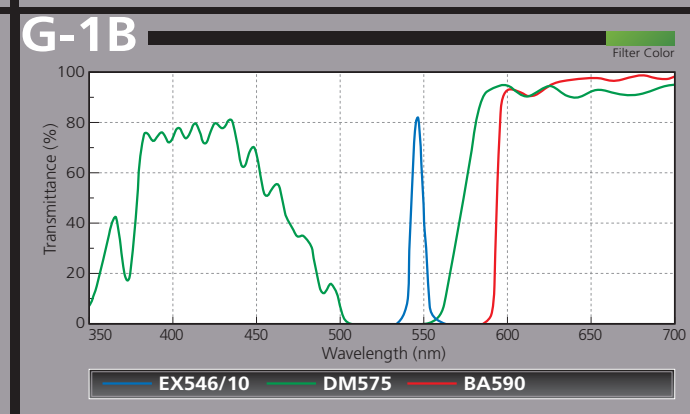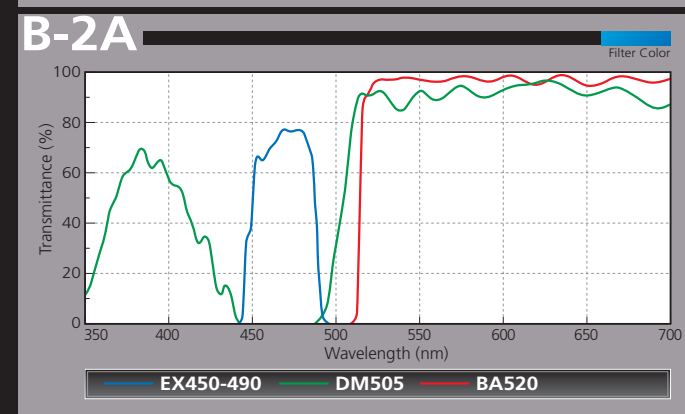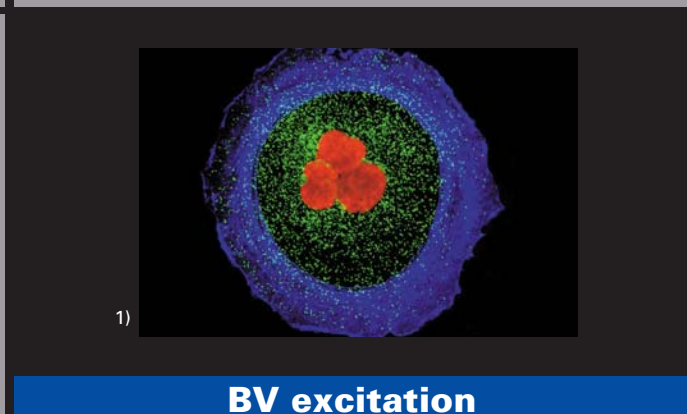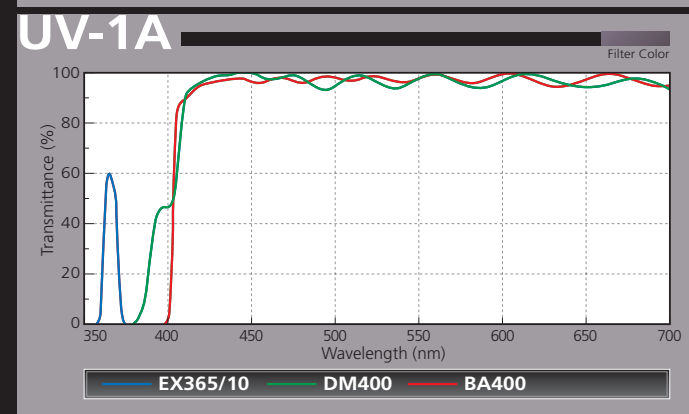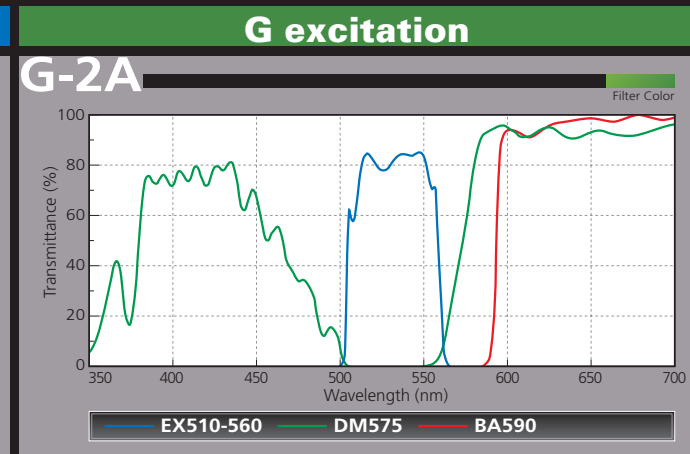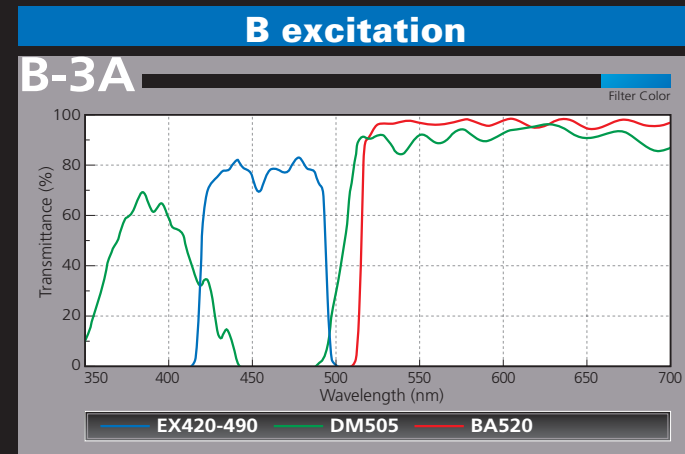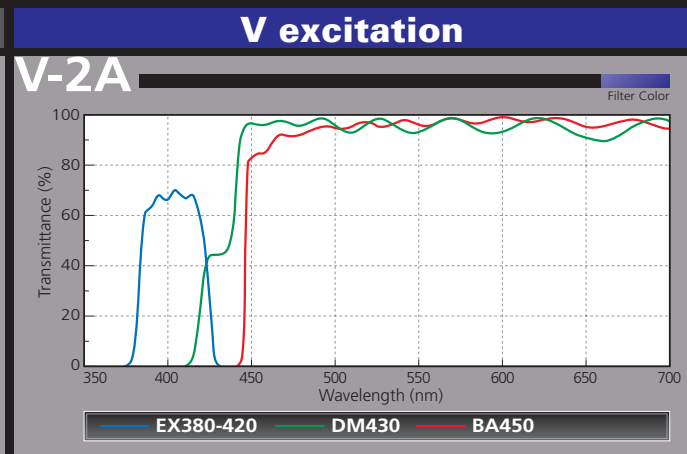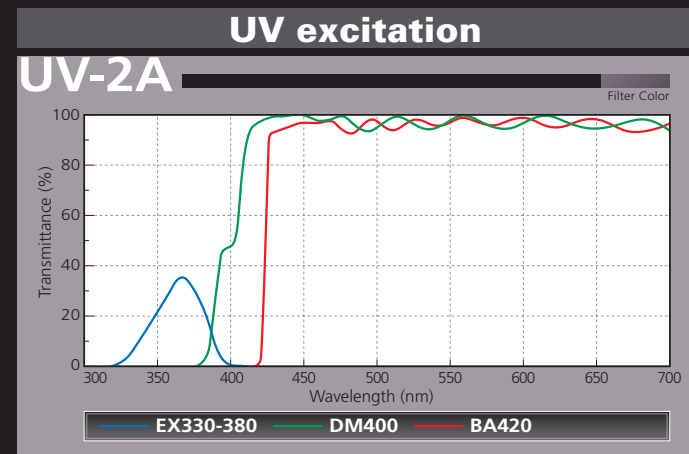

















EX: Excitation filter DM: Dichroic mirror BA: Barrier filter



# Fluorescence Filter Cubes from Nikon



## Special filter cubes for fluorescent reagents and fluorescent proteins

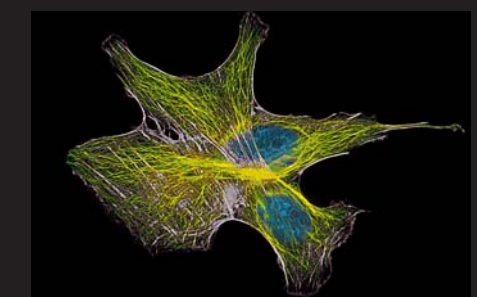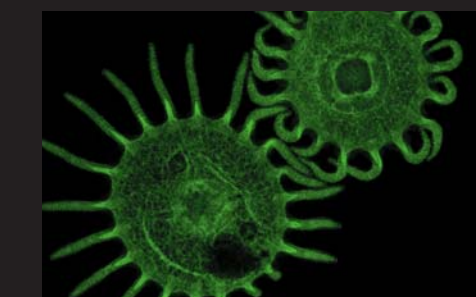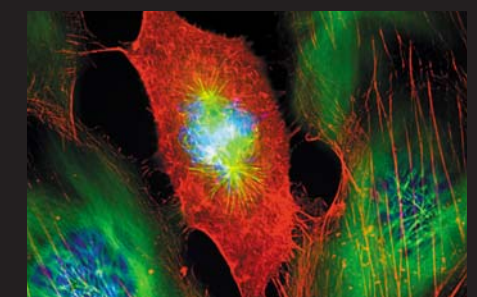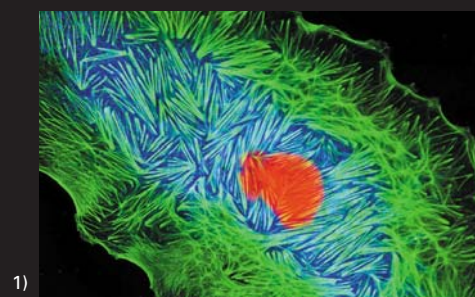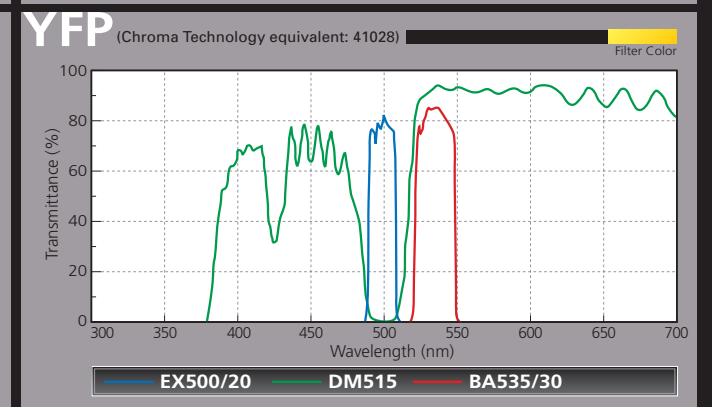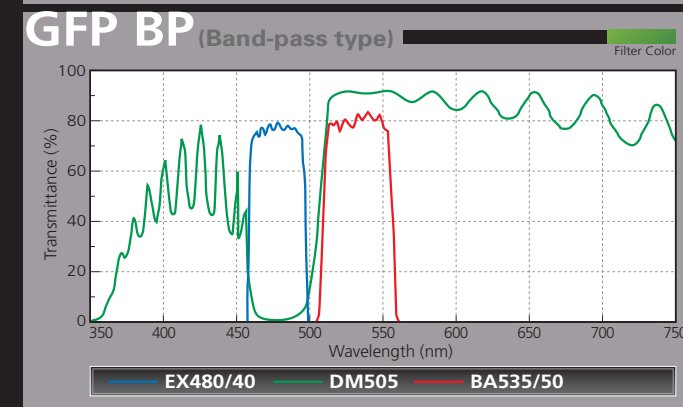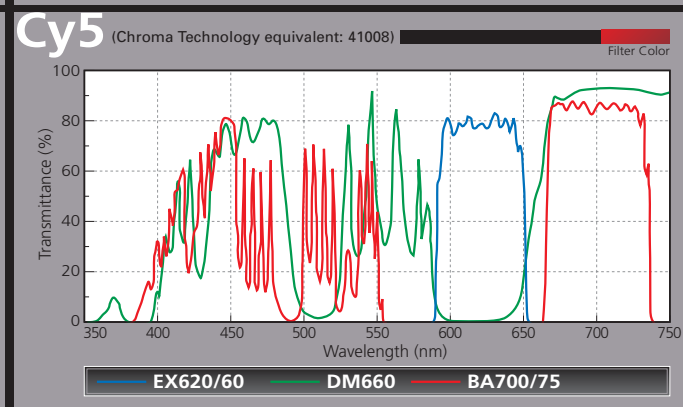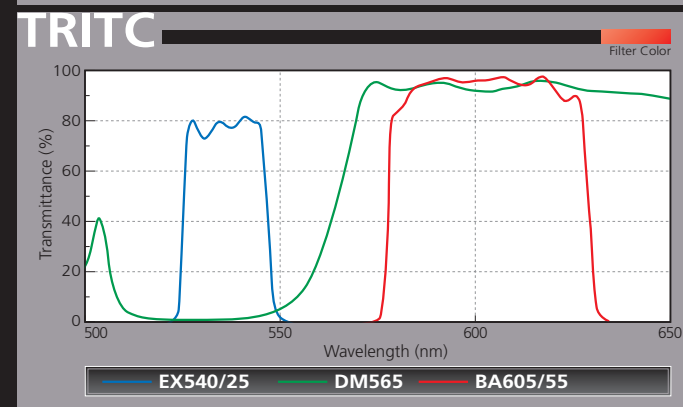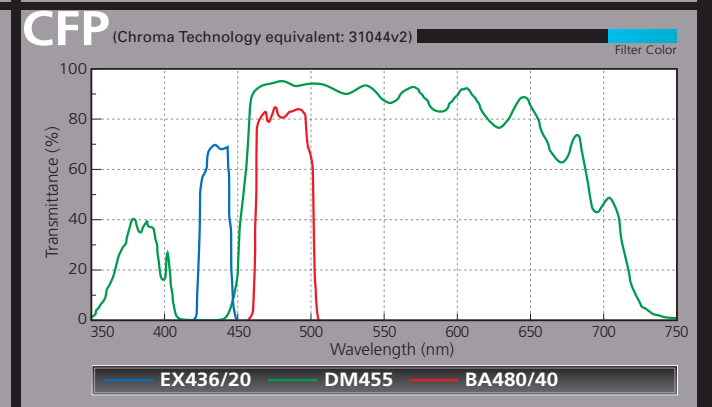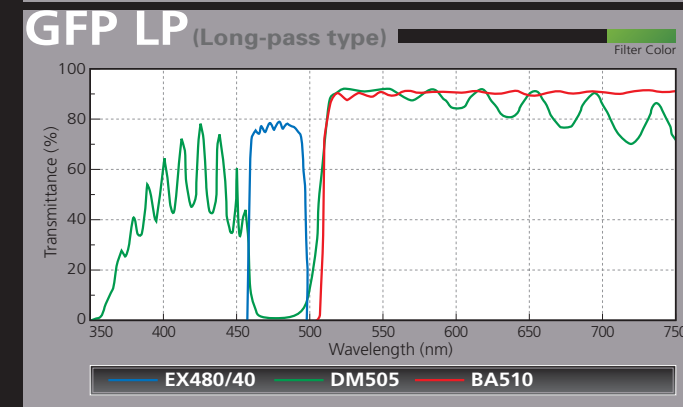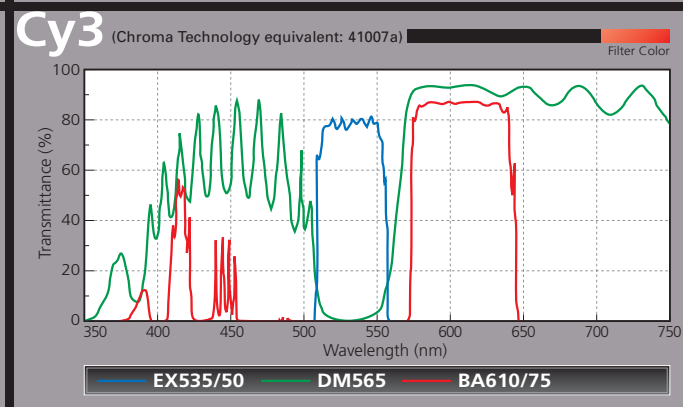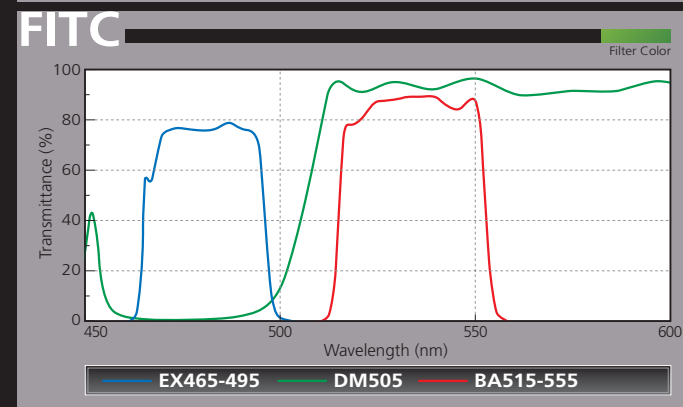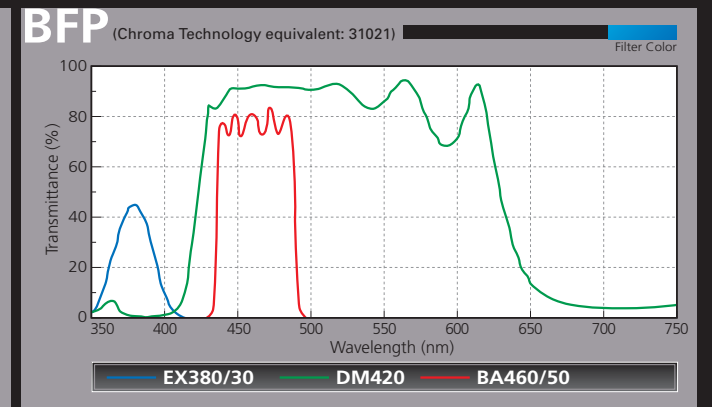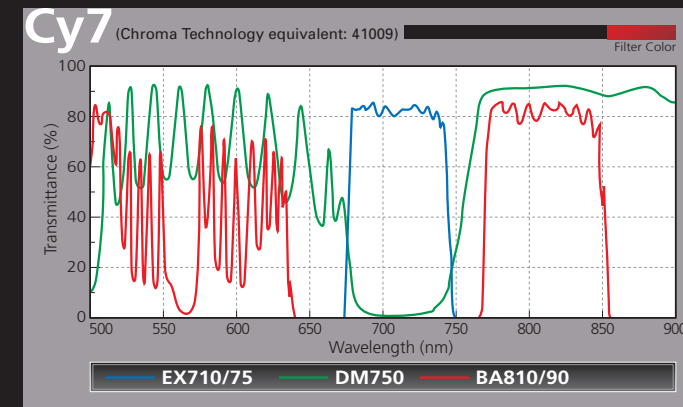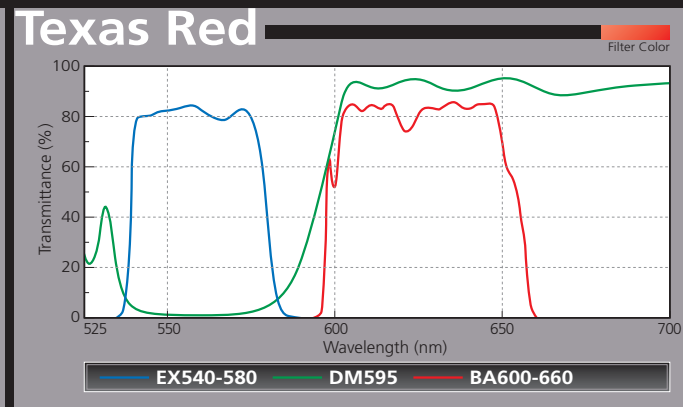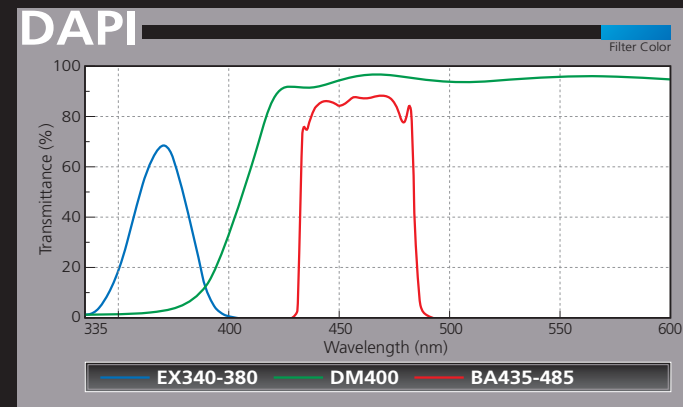







# Fluorescence Filter Cubes from Nikon



# Reagent Compatibility Table/ Specific Energy Distribution




## Multiband filter cubes

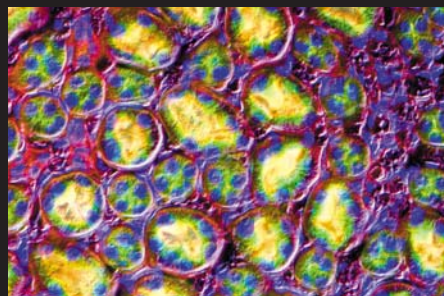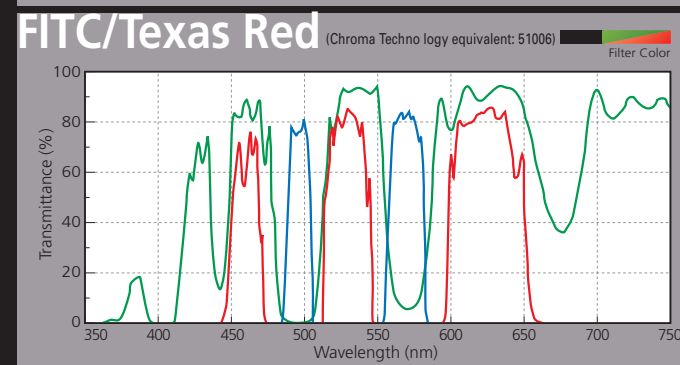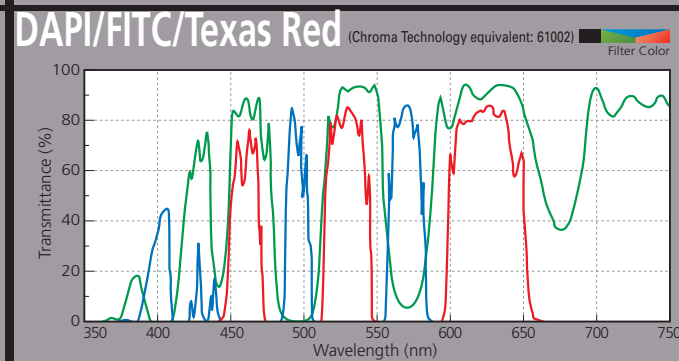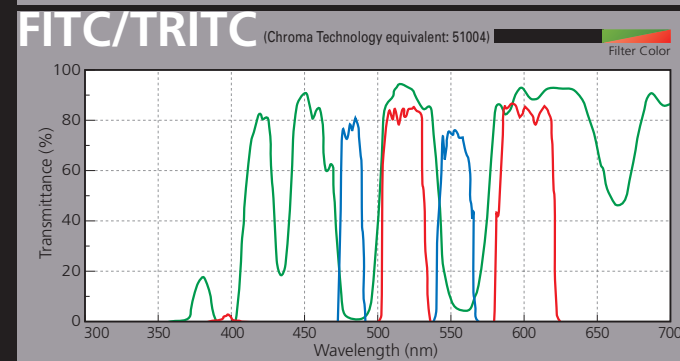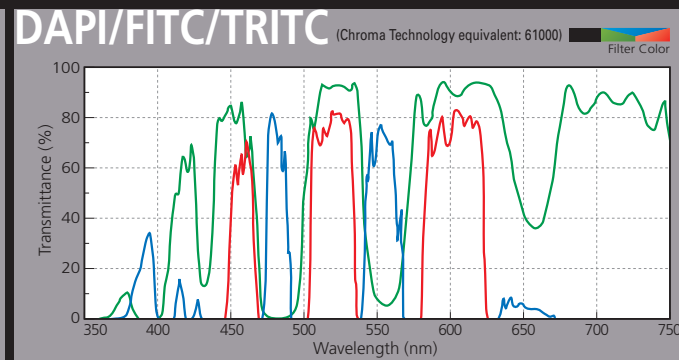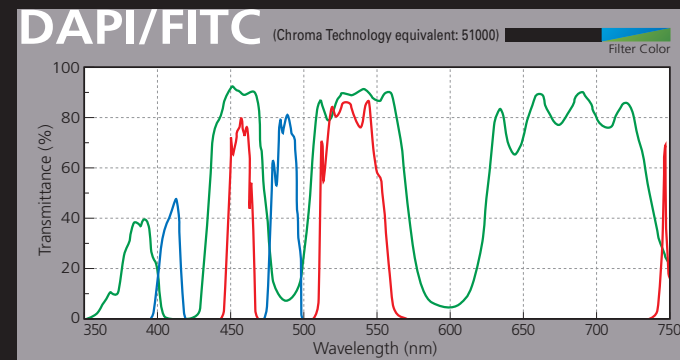

## Single filters can be switched for specific purposes

The optimal combination for the purpose of your observation can be created by easily removing the excitation filter, barrier filter, and/or the dichroic mirror.

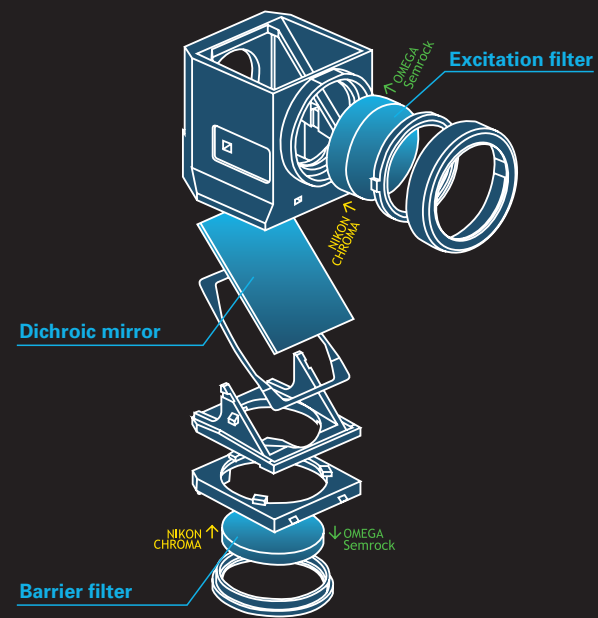


**Nikon and Chroma Technology:**  
The direction of the arrow on the excitation and barrier filter is the dichroic mirror side  
**Omega Optical and Semrock:**  
The direction of the arrow on the excitation/barrier filter is the direction of light

## Reagent compatibility table


Typical reagents	EX	EM	Nikon-compatible filter cubes
ACMA	430	474	BV-2B
Acridine Orange (DNA+RNA)	440-480	520-560	B-2A
Alexa Fluor 350	347	442	UV-2A
Alexa Fluor 488	495	519	B-2A, B-1A
Alexa Fluor 568	579	604	G-2A
Alexa Fluor 647	653	669	Cy5
Allophycocyanin (APC)	650	660	Cy5
BCECF (high ph)	503	528	B-2A
BFP (Blue Fluorescent Protein)	381	445	BFP
Calcein	494	517	FITC, B-2A
Calcium Green-1	506	531	B-2A
Cascade Blue	376	425	UV-2A
CFDA (Carboxyfluorescein)	495	520	FITC, B-2A
CFP (Cyan Fluorescent Protein)	458	480	CFP
Cy2	489	506	B-2A, GFP-BP
Cy3	550	570	Cy3, G-2A
Cy5	649	670	Cy5
DAPI	358	461	DAPI, UV-2E/C
DiOC6	480	501	B-2A, FITC-HYQ
Dil	549	565	Cy3, G-2A
DsRed (Red Fluorescent Protein)	558	583	TRITC, G-2E/C
Ethidium bromide	545	605	G-2A
FITC	494	518	FITC, B-2E/C
Fluo3	506	526	B-2A
FluoroGold	368	565	UV-2A
FM1-43	502	625	B-2A
Fura2	335	505	Fura-2
Fura Red	472	646	B-2A
Hoechst 33342 & 33258	352	461	UV-2A
Indo1	330	401	Indo-1
JC-1	514	529	B-2A, YFP
Lissamine rhodamine B	570	590	Cy3, G-2A, G-2B
Lusifer Yellow	428	536	B-3A
Lyso Tracker Green	505	511	B
MitoTracker Green	490	516	B-2A, FITC
MitoTracker Orange	551	576	Cy3, G-2A
Monochlorobimane	380	461	UV-2A
NBD (amine)	460	534	FITC, B-2A
Nile Red	549	628	G
Pacific Blue	405	455	V-2A
R-phycoerythrin	480/546/565	578	Cy3, G-2A*B-2A
POPO-3	534	570	Cy3, G-2A
Propidium iodide (PI)	536	617	G-2A
Pyronine Y	555	580	Cy3, G-2A
RH795	530	712	Cy3, G-2A
Rhodamine123	507	529	B-2A, FITC
SYTOX	504	523	B-2A, FITC
Texas Red	577	620	Texas Red, Y-2E/C
TMR (Tetramethylrhodamine)	555	580	TRITC, G-2E/C, Cy3
TO-PRO-3	642	661	Cy5
TOTO-3	642	660	Cy5
XRITC (X-rhodamine-5)	580	605	Texas Red, Cy3, Y-2E/C
YFP (Yellow Fluorescent Protein)	513	527	YFP
YOYO-1	491	509	B-2A, FITC, GFP-BP

## Specific energy distribution

### Mercury lamp

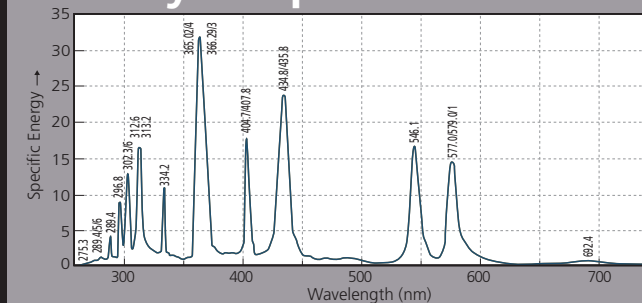

### Xenon lamp

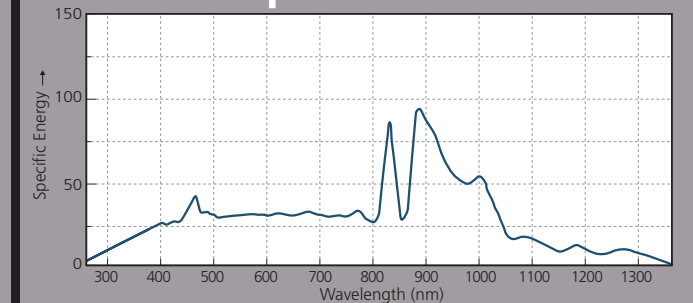






Qdot® conjugates have several special features, including extremely slow color fading, and they are winning acclaim as a new labeling tool for fluorescence observation. Nikon now offers a dedicated Qdot® detection filter, which maximizes the performance of the fluorescent probe.

## Qdot® nanocrystals

The quantum dot conjugate is made from nanometer-scale crystals of semiconductor material, and the color of light that they emit differs depending on the particle size. Qdot® conjugates are nanocrystals for labeling biomolecules such as antibodies and streptavidin. Unlike conventional organic dyes, Qdot® offers the following advantages:

- (1) Extremely slow color fading and long-term photo stability
- (2) Extremely high fluorescence intensity
- (3) Extremely sharp detection of wavelength distribution, enabling the simultaneous detection of different colors without any overlap
- (4) Compatible with all manner of optical microscopes, including confocal models


Absorption profiles (solid line) and emission profiles (dotted line) of Qdot® nanocrystals

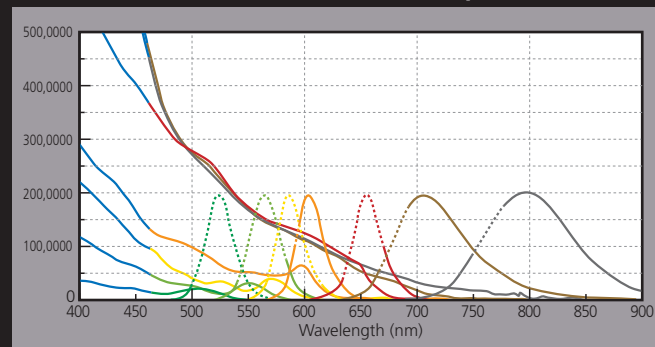

Qdot®525 conjugate excitation and emission	Green
Qdot®565 conjugate excitation and emission	Chartreuse
Qdot®585 conjugate excitation and emission	Yellow
Qdot®605 conjugate excitation and emission	Orange
Qdot®655 conjugate excitation and emission	Dark red
Qdot®705 conjugate excitation and emission	Near IR
Qdot®800(single/multi) conjugate excitation and emission	Near IR

Note: The emission of Qdot® 705 and 800 is not visible to the naked eye and must be detected with an IR-sensitive detector.

## Multi-color (5-color) immunostaining of fixed human epithelial cells

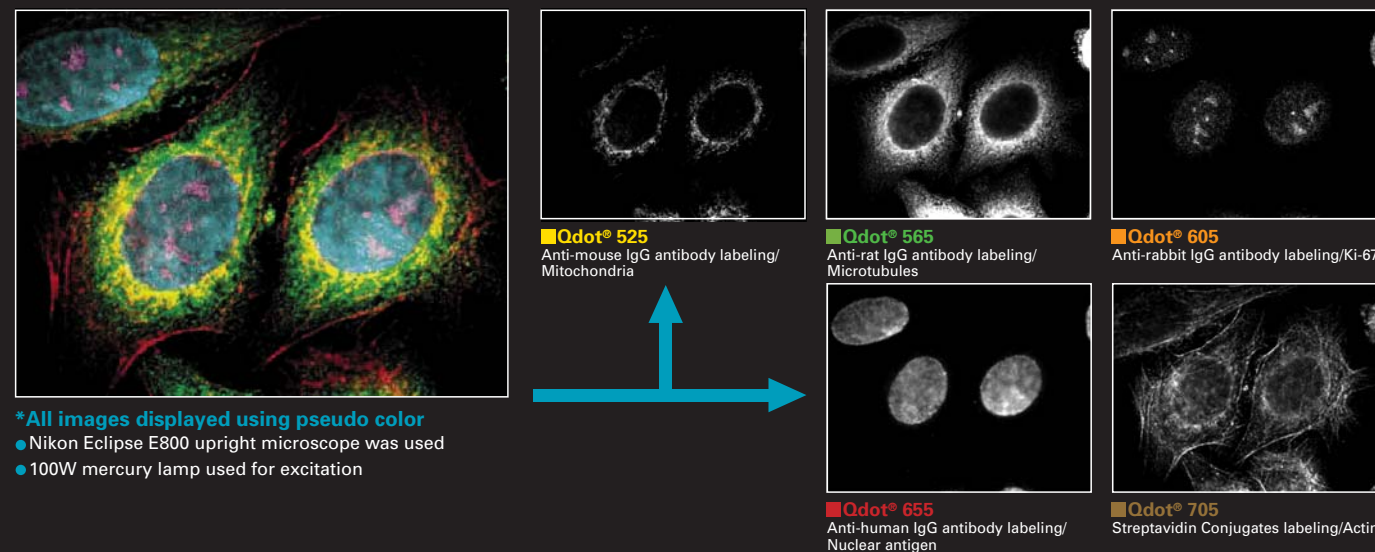

- \*All images displayed using pseudo color
- Nikon Eclipse E800 upright microscope was used
- 100W mercury lamp used for excitation


\* For Qdot® products, we recommend that the excitation area be changed according to your application.  
 Exciter 1: For researchers observing living cells (excitation at 402.5 nm to 447.5 nm) => D420/40  
 Exciter 2: For researchers observing fixed cells or who emphasize brightness and are not concerned about UV toxicity (excitation at 365 nm to 465 nm) => E460SPUV  
 Nikon also carries Qdot®-compatible filters for the C1 plus confocal laser microscope.

Select fluorescence filters and mirrors according to the following procedures.

## Checking the wavelength properties of fluorescent substances

Feel free to refer to the characteristic listed in catalogs and other sources to check the wavelength properties of fluorescent substances, but since these properties change slightly depending upon solution conditions, use a fluorometer or similar instrument to compare excitation and fluorescence wavelength properties while selecting the optimal optical element.

## Selecting a dichroic mirror

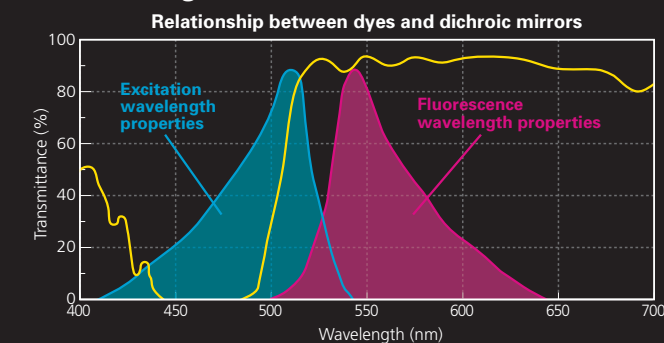

When comparing the spectral property curve of a dichroic mirror measured at a 45° angle, select a mirror for which the rising wavelength in the transmission range is in between the excitation wavelength and fluorescent wavelength of the fluorescent substance. When the excitation wavelength and fluorescent wavelength of the fluorescent substance are close to each other, select a dichroic mirror that rises closer to a short wavelength in order to transmit the fluorescence signal as much as possible. With a dichroic mirror, it is necessary to pay attention not only to the rising wavelength, but also the inclination of the rise. A dichroic mirror with a gentle rising edge may end up lowering the transmittance of the fluorescence signal and produce an unwanted crossover fluorescence emission signal.

## Selecting an barrier filter

Generally, a long-pass filter that transmits up to long wavelengths will be selected, but when observing a multi-stained specimen using discrete wavelength filter cubes, or when using a camera with high sensitivity in the high red or IR wavelength bands, select a band-pass filter that does not transmit the longer wavelengths.

## Selecting an excitation filter

Select a filter that satisfies the excitation wavelength of the fluorescent substance. Especially when using a mercury lamp as a light source, making the most of the emission lines of the lamp into the excitation wavelength enables highly efficient excitation. Since applying intense light in a wavelength range other than the excitation wavelength of the fluorescent substance being observed causes the background noise to rise and unnecessarily damages the sample, this situation should be avoided whenever possible.

## Combining an excitation and an barrier filter

When the wavelength properties of an excitation filter and barrier filter overlap, ideally, no light at all will pass through. Since fluorescence is a weak beam, even the slightest bit of light leakage will cause the background noise to rise, inviting degraded image quality.

Note: The wavelength characteristics of filters described in this catalogue are examples only. Filters may not fully match the specified wavelengths due to slight variations between production lots.